



#7241 Store at +4C

PathScan® Phospho-Met (Tyr1003) Sandwich ELISA Kit

Orders: 877-616-CELL (2355)
orders@cellsignal.com

Support: 877-678-TECH (8324)

Web: info@cellsignal.com
cellsignal.com

Species Cross Reactivity: H
UniProt ID: #P08581
Entrez-Gene Id: #4233

3 Trask Lane | Danvers | Massachusetts | 01923 | USA

For Research Use Only. Not for Use in Diagnostic Procedures.

Product Includes	Product #	Quantity	Color	Storage Temp
Phospho-Met (Tyr1003) Rabbit Detection mAb	13725	1 ea	Green (Lyophilized)	+4C
Anti-mouse IgG, HRP-linked Antibody (ELISA Formulated)	13304	1 ea	Red (Lyophilized)	+4C
Detection Antibody Diluent	13339	11 ml	Green	+4C
HRP Diluent	13515	11 ml	Red	+4C
TMB Substrate	7004	11 ml	Colorless	+4C
STOP Solution	7002	11 ml	Colorless	+4C
Sealing Tape	54503	2 ea		+4C
ELISA Wash Buffer (20X)	9801	25 ml	Colorless	+4C
ELISA Sample Diluent	11083	25 ml	Blue	+4C
Cell Lysis Buffer (10X)	9803	15 ml	Yellowish	-20C

Kit contents scale proportionally with size, except sealing tape.

Example: The V1 kit contains 5X the listed quantities above, but will exclude the sealing tape.

The microwell plate is supplied as 12 8-well modules - Each module is designed to break apart for 8 tests.

Description

The PathScan® Phospho-Met (Tyr1003) Sandwich ELISA Kit is a solid phase sandwich enzyme-linked immunosorbent assay (ELISA) that detects endogenous levels of Met when phosphorylated at Tyr1003. A Met Mouse Antibody has been coated onto the microwells. After incubation with cell lysates, Met (phospho and nonphospho) is captured by the coated antibody. Following extensive washing, a Phospho-Met (Tyr1003) Rabbit Detection Antibody is added to detect tyrosine phosphorylation of the captured Met protein. Anti-rabbit IgG, HRP-Linked Antibody is then used to recognize the bound detection antibody. HRP substrate, TMB, is added to develop color. The magnitude of the absorbance for this developed color is proportional to the quantity of Met phosphorylated at Tyr1003. Antibodies in kit are custom formulations specific to kit.

Specificity/Sensitivity

CST's PathScan® Phospho-Met (Tyr1003) Sandwich ELISA Kit #7241 detects Met when phosphorylated on Tyr1003. By using this kit, high levels of phospho-Met protein at Tyr1003 are detected in HCC827 cells where Met is constitutively phosphorylated (Figure 1). These high levels are abolished in HCC827 cells lysed without addition of phosphatase inhibitors* to the lysis buffer. The levels of total Met protein (phospho and nonphospho) as detected by PathScan® Total Met Sandwich ELISA Kit #7242 remain unchanged (Figure 1).

* Phosphatase inhibitors include sodium pyrophosphate, β -glycerophosphate and Na_3VO_4 .

This kit detects proteins from the indicated species, as determined through in-house testing, but may also detect homologous proteins from other species.

Background

Met, a high affinity tyrosine kinase receptor for hepatocyte growth factor (HGF, also known as scatter factor) is a disulfide-linked heterodimer made of 45 kDa α - and 145 kDa β -subunits (1,2). The α -subunit and the amino-terminal region of the β -subunit form the extracellular domain. The remainder of the β -chain spans the plasma membrane and contains a cytoplasmic region with tyrosine kinase activity. Interaction of Met with HGF results in autophosphorylation at multiple tyrosines, which recruit several downstream signaling components, including Gab1, c-Cbl, and PI3 kinase (3). These fundamental events are important for all of the biological functions involving Met kinase activity. The addition of a phosphate at cytoplasmic Tyr1003 is essential for Met protein ubiquitination and degradation (4). Phosphorylation at Tyr1234/1235 in the Met kinase domain is critical for kinase activation. Phosphorylation at Tyr1349 in the Met cytoplasmic domain provides a direct binding site for Gab1 (5). Research studies have shown that altered Met levels and/or tyrosine kinase activities are found in several types of tumors, including renal, colon, and breast. Thus, investigators have concluded that Met is an attractive potential cancer therapeutic and diagnostic target (6,7).

Background References

1. Cooper, C.S. et al. (1984) *Nature* 311, 29-33.
 2. Bottaro, D.P. et al. (1991) *Science* 251, 802-4.
 3. Bardelli, A. et al. (1997) *Oncogene* 15, 3103-11.
 4. Taher, T.E. et al. (2002) *J Immunol* 169, 3793-800.
 5. Schaeper, U. et al. (2000) *J Cell Biol* 149, 1419-32.
 6. Eder, J.P. et al. (2009) *Clin Cancer Res* 15, 2207-14.
 7. Sattler, M. and Salgia, R. (2009) *Update Cancer Ther* 3, 109-118.
-

Trademarks and Patents

Cell Signaling Technology is a trademark of Cell Signaling Technology, Inc.

PathScan is a registered trademark of Cell Signaling Technology, Inc.

All other trademarks are the property of their respective owners. Visit cellsignal.com/trademarks for more information.

Limited Uses

Except as otherwise expressly agreed in a writing signed by a legally authorized representative of CST, the following terms apply to Products provided by CST, its affiliates or its distributors. Any Customer's terms and conditions that are in addition to, or different from, those contained herein, unless separately accepted in writing by a legally authorized representative of CST, are rejected and are of no force or effect.

Products are labeled with For Research Use Only or a similar labeling statement and have not been approved, cleared, or licensed by the FDA or other regulatory foreign or domestic entity, for any purpose. Customer shall not use any Product for any diagnostic or therapeutic purpose, or otherwise in any manner that conflicts with its labeling statement. Products sold or licensed by CST are provided for Customer as the end-user and solely for research and development uses. Any use of Product for diagnostic, prophylactic or therapeutic purposes, or any purchase of Product for resale (alone or as a component) or other commercial purpose, requires a separate license from CST. Customer shall (a) not sell, license, loan, donate or otherwise transfer or make available any Product to any third party, whether alone or in combination with other materials, or use the Products to manufacture any commercial products, (b) not copy, modify, reverse engineer, decompile, disassemble or otherwise attempt to discover the underlying structure or technology of the Products, or use the Products for the purpose of developing any products or services that would compete with CST products or services, (c) not alter or remove from the Products any trademarks, trade names, logos, patent or copyright notices or markings, (d) use the Products solely in accordance with CST Product Terms of Sale and any applicable documentation, and (e) comply with any license, terms of service or similar agreement with respect to any third party products or services used by Customer in connection with the Products.

#7241

PathScan® Phospho-Met (Tyr1003) Sandwich ELISA Kit



ELISA Colorimetric

NOTE: Refer to product-specific datasheets or product webpage for assay incubation temperature.

A. Solutions and Reagents

NOTE: Prepare solutions with reverse osmosis deionized (RODI) or equivalent grade water.

1. **20X Phosphate Buffered Saline (PBS):** (#9808) To prepare 1 L PBS: add 50 ml 10X PBS to 950 ml dH₂O, mix.
2. Bring all microwell strips to room temperature before use.
3. Prepare 1X Wash Buffer by diluting 20X Wash Buffer (included in each PathScan® Sandwich ELISA Kit) in dH₂O.
4. **1X Cell Lysis Buffer:** 10X Cell Lysis Buffer (#9803): To prepare 10 ml of 1X Cell Lysis Buffer, add 1 ml of 10X Cell Lysis Buffer to 9 ml of dH₂O, mix. Buffer can be stored at 4°C for short-term use (1-2 weeks).

Recommended: Add 1 mM phenylmethylsulfonyl fluoride (PMSF) (#8553) immediately before use.

NOTE: Refer to product-specific datasheet or webpage for lysis buffer recommendation.

5. **TMB Substrate:** (#7004).
6. **STOP Solution:** (#7002).

B. Preparing Cell Lysates

For adherent cells

1. Aspirate media when the culture reaches 80–90% confluence. Treat cells by adding fresh media containing regulator for desired time.
2. Remove media and rinse cells once with ice-cold 1X PBS.
3. Remove PBS and add 0.5 ml ice-cold 1X cell lysis buffer plus 1 mM PMSF to each plate (10 cm diameter) and incubate the plate on ice for 5 min.
4. Scrape cells off the plate and transfer to an appropriate tube. Keep on ice.
5. Sonicate lysates on ice.
6. Microcentrifuge for 10 min (x14,000 rpm) at 4°C and transfer the supernatant to a new tube. The supernatant is the cell lysate. Store at -80°C in single-use aliquots.

For suspension cells

1. Remove media by low speed centrifugation (~1,200 rpm) when the culture reaches 0.5–1.0 x 10⁶ viable cells/ml. Treat cells by adding fresh media containing regulator for desired time.
2. Collect cells by low speed centrifugation (~1,200 rpm) and wash once with 5–10 ml ice-cold 1X PBS.
3. Cells harvested from 50 ml of growth media can be lysed in 2.0 ml of 1X cell lysis buffer plus 1 mM PMSF.
4. Sonicate lysates on ice.
5. Microcentrifuge for 10 min (x14,000 rpm) at 4°C and transfer the supernatant to a new tube. The supernatant is the cell lysate. Store at -80°C in single-use aliquots.

C. Test Procedure

1. After the microwell strips have reached room temperature, break off the required number of microwells. Place the microwells in the strip holder. Unused microwells must be resealed in the storage bag and stored at 4°C immediately.
2. Cell lysates can be undiluted or diluted with sample diluent (supplied in each PathScan® Sandwich ELISA Kit, blue color). Individual datasheets or product webpage for each kit provide information regarding an appropriate dilution factor for lysates and kit assay results.
3. Add 100 µl of each undiluted or diluted cell lysate to the appropriate well. Seal with tape and press firmly onto top of microwells. Incubate the plate for 2 hr at 37°C. Alternatively, the plate can be incubated overnight at 4°C.
4. **Gently remove the tape and wash wells:**
 1. Discard plate contents into a receptacle.
 2. Wash 4 times with 1X wash buffer, 200 µl each time per well.
 3. For each wash, strike plates on fresh paper towels hard enough to remove the residual solution in each well, but do not allow wells to completely dry at any time.
 4. Clean the underside of all wells with a lint-free tissue.
5. Add 100 µl of detection antibody (green color) to each well. Seal with tape and incubate the plate at 37°C for 1 hr.

6. Repeat wash procedure (Section C, Step 4).
7. Add 100 μ l of HRP-linked secondary antibody (red color) to each well. Seal with tape and incubate the plate for 30 min at 37°C.
8. Repeat wash procedure (Section C, Step 4).
9. Add 100 μ l of TMB substrate to each well. Seal with tape and incubate the plate for 10 min at 37°C or 30 min at 25°C.
10. Add 100 μ l of STOP solution to each well. Shake gently for a few seconds.

NOTE: Initial color of positive reaction is blue, which changes to yellow upon addition of STOP solution.

11. Read results

1. **Visual Determination:** Read within 30 min after adding STOP solution.
2. **Spectrophotometric Determination:** Wipe underside of wells with a lint-free tissue. Read absorbance at 450 nm within 30 min after adding STOP solution.

posted June 2005

revised November 2013