

Vimentin Antibody Sampler Kit



Orders: 877-616-CELL (2355)
orders@cellsignal.com

Support: 877-678-TECH (8324)

Web: info@cellsignal.com
cellsignal.com

3 Trask Lane | Danvers | Massachusetts | 01923 | USA

1 Kit (3 x 20 microliters)

For Research Use Only. Not for Use in Diagnostic Procedures.

Product Includes	Product #	Quantity	Mol. Wt	Isotype/Source
Phospho-Vimentin (Ser56) Antibody	3877	20 µl	57 kDa	Rabbit
Phospho-Vimentin (Ser83) Antibody	3878	20 µl	57 kDa	Rabbit
Vimentin (D21H3) XP [®] Rabbit mAb	5741	20 µl	57 kDa	Rabbit IgG
Anti-rabbit IgG, HRP-linked Antibody	7074	100 µl		Goat

Please visit cellsignal.com for individual component applications, species cross-reactivity, dilutions, protocols, and additional product information.

Description

The Vimentin Antibody Sampler Kit provides an economical means to detect total levels of vimentin, vimentin phosphorylated at Ser56, and vimentin phosphorylated at Ser83. The kit contains enough primary and secondary antibody to perform two western blot experiments.

Storage

Supplied in 10 mM sodium HEPES (pH 7.5), 150 mM NaCl, 100 µg/ml BSA, 50% glycerol and less than 0.02% sodium azide. Store at -20°C. Do not aliquot the antibody.

Background

The cytoskeleton consists of three types of cytosolic fibers: microfilaments (actin filaments), intermediate filaments, and microtubules. Major types of intermediate filaments are distinguished by their cell-specific expression: cytokeratins (epithelial cells), glial fibrillary acidic protein (GFAP) (glial cells), desmin (skeletal, visceral, and certain vascular smooth muscle cells), vimentin (mesenchyme origin), and neurofilaments (neurons). GFAP and vimentin form intermediate filaments in astroglial cells and modulate their motility and shape (1). In particular, vimentin filaments are present at early developmental stages, while GFAP filaments are characteristic of differentiated and mature brain astrocytes. Thus, GFAP is commonly used as a marker for intracranial and intraspinal tumors arising from astrocytes (2). Research studies have shown that vimentin is present in sarcomas, but not carcinomas, and its expression is examined in conjunction with that of other markers to distinguish between the two (3). Vimentin's dynamic structural changes and spatial re-organization in response to extracellular stimuli help to coordinate various signaling pathways (4). Phosphorylation of vimentin at Ser56 in smooth muscle cells regulates the structural arrangement of vimentin filaments in response to serotonin (5,6). Remodeling of vimentin and other intermediate filaments is important during lymphocyte adhesion and migration through the endothelium (7). During mitosis, CDK1 phosphorylates vimentin at Ser56. This phosphorylation provides a PLK binding site for vimentin-PLK interaction. PLK further phosphorylates vimentin at Ser83, which might serve as memory phosphorylation site and play a regulatory role in vimentin filament disassembly (8,9). Additionally, studies using various soft-tissue sarcoma cells have shown that phosphorylation of vimentin at Ser39 by Akt1 enhances cell migration and survival, suggesting that vimentin could be a potential target for soft-tissue sarcoma targeted therapy (10,11).

Background References

- Eng, L.F. et al. (2000) *Neurochem Res* 25, 1439-51.
- Goebel, H.H. et al. (1987) *Acta Histochem Suppl* 34, 81-93.
- Leader, M. et al. (1987) *Histopathology* 11, 63-72.
- Helfand, B.T. et al. (2004) *J Cell Sci* 117, 133-41.
- Tang, D.D. et al. (2005) *Biochem J* 388, 773-83.
- Fomina, I.G. et al. (1990) *Klin Med (Mosk)* 68, 125-7.
- Nieminen, M. et al. (2006) *Nat Cell Biol* 8, 156-62.
- Yamaguchi, T. et al. (2005) *J Cell Biol* 171, 431-6.
- Oguri, T. et al. (2006) *Genes Cells* 11, 531-40.
- Zhu, Q.S. et al. (2011) *Oncogene* 30, 457-70.
- Xue, G. and Hemmings, B.A. (2013) *J Natl Cancer Inst* 105, 393-404.

Trademarks and Patents

Cell Signaling Technology is a trademark of Cell Signaling Technology, Inc.

All other trademarks are the property of their respective owners. Visit cellsignal.com/trademarks for more information.

Limited Uses

Except as otherwise expressly agreed in a writing signed by a legally authorized representative of CST, the following terms apply to Products provided by CST, its affiliates or its distributors. Any Customer's terms and conditions that are in addition to, or different from, those contained herein, unless

separately accepted in writing by a legally authorized representative of CST, are rejected and are of no force or effect.

Products are labeled with For Research Use Only or a similar labeling statement and have not been approved, cleared, or licensed by the FDA or other regulatory foreign or domestic entity, for any purpose. Customer shall not use any Product for any diagnostic or therapeutic purpose, or otherwise in any manner that conflicts with its labeling statement. Products sold or licensed by CST are provided for Customer as the end-user and solely for research and development uses. Any use of Product for diagnostic, prophylactic or therapeutic purposes, or any purchase of Product for resale (alone or as a component) or other commercial purpose, requires a separate license from CST. Customer shall (a) not sell, license, loan, donate or otherwise transfer or make available any Product to any third party, whether alone or in combination with other materials, or use the Products to manufacture any commercial products, (b) not copy, modify, reverse engineer, decompile, disassemble or otherwise attempt to discover the underlying structure or technology of the Products, or use the Products for the purpose of developing any products or services that would compete with CST products or services, (c) not alter or remove from the Products any trademarks, trade names, logos, patent or copyright notices or markings, (d) use the Products solely in accordance with CST Product Terms of Sale and any applicable documentation, and (e) comply with any license, terms of service or similar agreement with respect to any third party products or services used by Customer in connection with the Products.

Orders: 877-616-CELL (2355) • orders@cellsignal.com • Support: 877-678-TECH (8324) • info@cellsignal.com • Web: cellsignal.com
For Research Use Only. Not for Use in Diagnostic Procedures.